
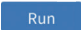


ANCHOR

SpeedPURE® Flex Kit (384) (RUO)

- On the display select  (upper left corner)
- Choose the desired protocol (Fig. 4)
 - SpeedPURE_200 for 200 µL sample input
 - SpeedPURE_400 for 400 µL sample input
- Select  (lower right corner) to start the sample preparation
- After the run is complete:
 - Remove and discard magnetic rod covers
 - Remove Deep Well Plates from instrument and collect eluates from columns 6 and 12 (Fig. 5)
 - Discard used Deep Well Plates

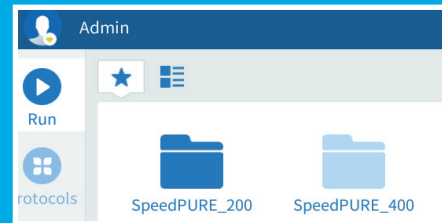
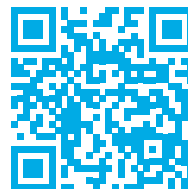


Fig. 4

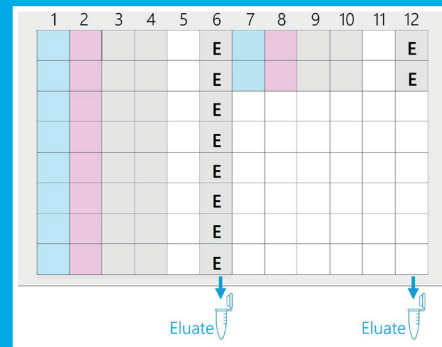


Fig. 5

REF AP03400L **QG** AP03491L-EN, 2026-02-21

Contact

ANCHOR Diagnostics GmbH
Grandweg 64 | D-22529 Hamburg

Support

+49 40 52 471 62 0
support@anchor-diagnostics.com